## Meßmikroskop Measuring microscope *Microscope de mesure*

### MM-KKE-M-C-U

Typenschlüssel / Model code / *Code de commande*

MM KKE M C U

**Grundtyp** / Basic model / *Modèle de base*

MM = **Meßmikroskop** / Measuring microscope / *Microscope de mesure*

**Optik** / Lens system / *Optique*

KKE = **Triokular** / Triocular / *Triculaire*

**Versorgungsspannung** / Supply voltage / *Tension d'alimentation*

0 = 240 V 50 Hz (Australien)

M = 230 V 50 Hz (Europa)

P = 110 V 60 Hz (Japan)

**Zubehör** / Accessories / *Accessoires*

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**Bilddigitalisierung** / Image digitization / *Numérisation de l'image*

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CCD camera with USB port to laptop or PC

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### Product description

This measuring microscope is mainly designed to measure particles on filter membranes of oil samples.

The microscope is sturdy and stable in its construction and offers a high level of operating comfort. The body is adjusted by an easily operated tangent screw for coarse adjustment and a fine adjustment tangent screw, thus ensuring optimum focus at maximum magnification. The integrated lighting with a plug-in power supply unit provides sufficient image brightness even with 200x magnification. The stand is fitted with a 3-part knurled revolving unit and a slide table with X and Y movements. The optical fittings consist of the achromatic lenses: 4:1, 10:1, 20:1. These lenses are used in combination with a micrometer eyepiece with 10x magnification. The measuring scale in the eyepiece has a scale division of 1 mm in 100 parts. By using the micrometer eyepiece and the enclosed measuring card, it is possible to determine the size of the subject directly with all three lenses. The microscope cabinet protects the microscope from knocks and dust.

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## NOTE

The information in this brochure relates to the operating conditions and applications described.

For applications or operating conditions not described, please contact the relevant technical department.

Subject to technical modifications.

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